



U ■ ○ ■ //

U

—

—

—

—



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
Copyright © by Bruker Corporation

All rights reserved. No part of this publication may be reproduced, stored in a retrieval system, or transmitted, in any form, or by any means without the prior consent of the publisher. Product names used are trademarks or registered trademarks of their respective holders.

© September 08, 2020 Bruker Corporation

Document Number:

P/N: H179023



#	1	1	
%	1.1	How to Use this Manual	5
)			
+			
/	4.1	TopSpin Interface - Overview	7
	4.2	First Start of TopSpin - Load a Dataset	8
	4.3	Basic Processing and Manipulations	10
	4.4	Basic Processing	11
	4.5	Phase Correction	12
	4.6	Baseline Correction.....	13
	4.7	Axis Calibration	13
	4.8	Advanced Processing Options.....	13
(5.1	Peak Picking	14
	5.2	Spectrum Integration.....	15
	5.3	Further Analysis Options.....	16
.	6.1	Print Screen Directly	17
	6.2	Print Spectrum Using Layout (Plot).....	18
1			
8			

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


U . O . /
~ i"/fil° k t Ž k a ž i"/"


U . O . /
~ i"/fil° k t ž k a ž i"/"


U . o . /
~ i " / f i n ° k t ž e n ž i " ! "






17. H1\$*~"1#~Z" /L~

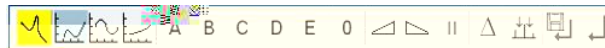For further improvement of the spectrum baseline,  **Baseline** leads to the interactive baseline correction. The dataset window will show icons specific for baseline correction.



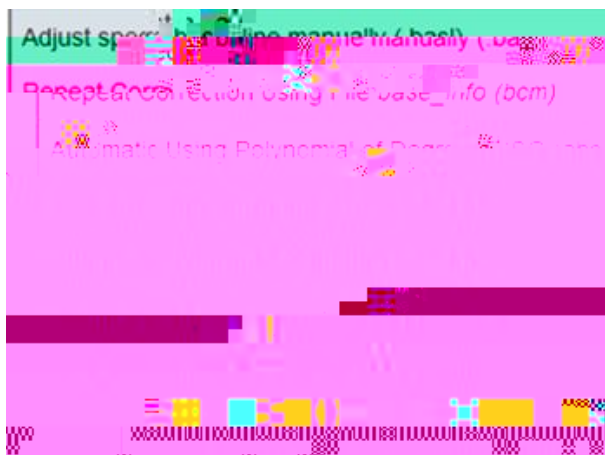
In addition, the Baseline pulldown menu gives access to further baseline correction operations.



17.1 21\$1#1*~JZL~



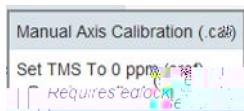
gives access to following pulldown menu with axis calibration options:




17.8 2!E1 /" !16Z~/"\$\$~ 01D~L~ \$



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gives access to further options through following pulldown menu:




(2 Ł*3\$ "ŹŁ4"fi° ' " /ŁŽ i 5

To enter the Analyse Process, click , the following workflow appears:



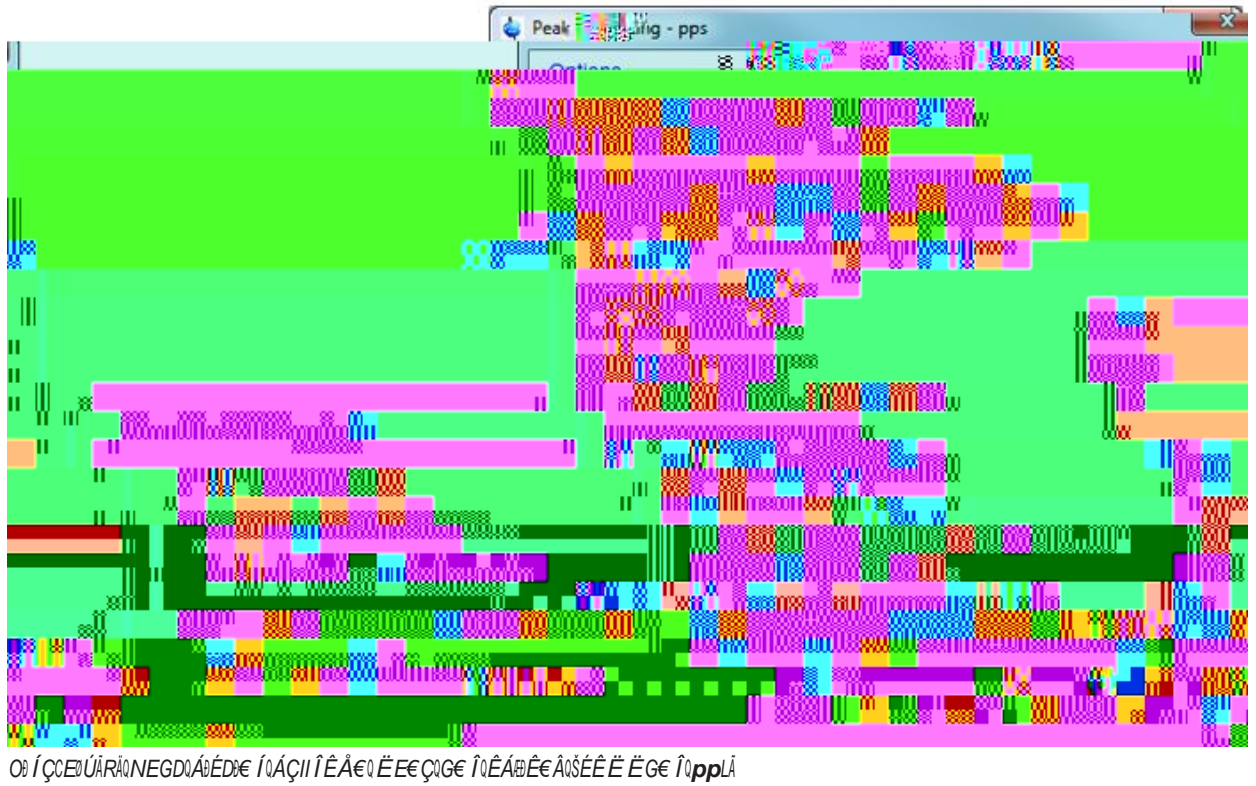
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
08 Í ÇCE0ÚÁPÁ0 Ú ÊCDFIÊÁ0BÇÊÊ€0BGCQFÊC0ÁÁEÊ€ÇÊ0G€GIYÁ0ÁÁ

(*% 6" Łfil6"/fi" 0

 gives access to the interactive peak picking. The dataset window shows specific icons for peak picking:




In addition, the Pick Peaks pulldown menu gives access to further peak picking operations and options:



0i ÍÇEÜÁRÁ00€E ÍÇG0É€QÁÇII ÎÊÁ€0 ÊE€ÇQÇE ÍQÊÁBÊEÁ0SÊÊÊ ÊE€ ÎQpplÄ

(') ° · "/Łž i 5 fl& Ł "0žŁŁž ~

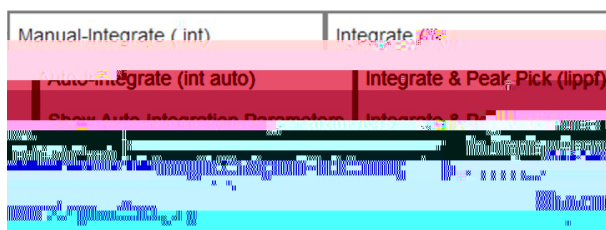
Integrate button gives access to the integration pulldown menu:



0i ÍÇEÜÁRÁ00€E ÍÇG0É€QÁÇII ÎÊÁ€0 ÊE€ÇQÇE ÍQ00€E ÍÇG0É€QÉ ÍQÊEÁQ00iE0 ÊE€ÇQŠÊÊÊ ÊE€ ÎQintrngLÄ

In addition, more options can be obtained through the command ~ Ł:

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
U . O . /
" i " / f i n ° k t ž e n ž i " ! "




• **6Ž" ĸ" 0fl4 "fl ° " /ŁŽ 1 5**


Printing the spectrum can be done in two different ways:

• '% **6Ž" ĸfl ° /Ž" " fl7Ž" /Ł*3**

Click on the printer icon  , located on the upper right part of the TopSpin window, allows direct printing of the spectrum. Details can be obtained as usual by dragging the mouse on the icon. Clicking the button gives access to the print window:

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

1) 6Z~_Łfl° · "/ŁŽ1 5 fl; \$~_0fl- Ł3~ 1Łfl>6*~Ł?

To use this TopSpin feature, click the button , or use the *~Ł command in the TopSpin command line. The following TopSpin window appears:



OB ÍÇÇEBŪĀRĀQNGCERĒPQ ĪĒĀ Ī ĀPĒQĀĪĒRQ.ĀPĒ ĪĒĀQŠ.ĀPĒQĀIGYĒÇELĀ

The left part of the plot window shows the names of the selected layout and printer. They can be changed with the arrow button, as well as other plotting view parameters, like limits or display. New NMR elements (title, parameters, 2D spectrum, etc.) can be inserted.

The right part presents the spectrum like it will be plotted. Changes on the left part will be shown in real time in the layout on the right.

1 , 1ŽŁ4"Žfl7~ / 1 5 " ŁŁŁ~

For detailed information please refer to the TopSpin Manual section under  =fl < Ł ĩ Ł* \$fl> ! ~ / \$?



8 # ~ ŁŁ/Ł

<Ł ŁŁŁ/ŁŁŹ"Ź

Bruker BioSpin GmbH
Silberstreifen 4
D-76287 Rheinstetten
Germany

E-Mail: € ĘHÂÇÁÁĘŒ Z BCÇDEÇĀĘĘĘ
AĘĘĀĀĀĀĀĀBCÇDEÇĀĘĘĘ
WEEE DE43181702

HŹ i fi"ŹfiH~°~ŁŁfi9~Ł*Ł"\$

Contact our Bruker BioSpin service centers.

HŽi fi"Ži#~Ž'~Ži~

info@bruker.com
www.bruker.com